



Optimice® unpacking instructions:

Copies of these instructions are also available for download in a PDF format at: www.animalcaresystems.com
 Pallet is accessible from two sides with a forklift or pallet jack. Be sure your equipment can handle the weight and size. USE CAUTION WHEN LIFTING RACKS!! RACKS ARE TOP HEAVY!! Rack weights and dimensions are noted below.
IF YOUR RACKS ARE SHIPPED VIA OCEAN CONTAINER (NOT BOXED) PLEASE SKIP TO STEP #5

TOOLS REQUIRED: Phillips driver, strap cutters and 7/16" (11mm) wrench.

1. Cut the vertical and horizontal straps.
2. Remove the cardboard top.
3. Remove the two large cardboard sides.
4. Remove the 4 cardboard/foam corners from the rack.

CAUTION: Check the corner pieces for any accessories that may be packed inside! They will be marked with yellow CAUTION TAPE.




5. Unscrew the two 1x6s from both sides of the pallet with a Phillips driver. Use these l



OptiMICE® Racks		
	Rack Only	Rack/pallet & box
	Lbs (Kg)	Lbs (Kg)
Standard Rack	560 (254)	590 (268)
Rack w/auto water	586 (266)	616 (279)
	Inches (mm)	Inches (mm)
Length	40 (1016)	44 (1118)
Width	34 (864)	44 (1118)
Height	78 (1981)	79 (2006)

6. Remove 4 hex bolts with 7/16" (or 11mm) wrench from underneath the unit at each corner near the casters.

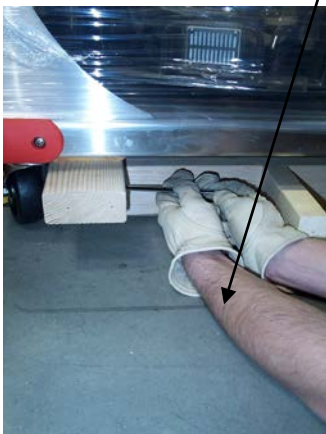

1 x 6 used as a pry bar

7. Using the 1x6 from above, pry up one side of the rack and slide out a 1x4. Then pry up the other side and remove the other 1x4.
8. OptiMICE® should now be on its wheels ready to roll off the remaining pallet.
9. Carefully remove the plastic wrap.