

Optimice SOP - Using Standard Cage Components

Feeder

- 1. Fill the feeder, ensuring the feed does not go beyond the upper rim. Otherwise, the cage top may not securely clip in place.
- 2. Place the feeder upon the hanger within the cage.





Water Bottle

- 1. After filling, cap the bottle; twist the cap one-quarter turn to secure the cap and prevent leaks.
- 2. Hang the bottle on the hook beside the feeder.
- 3. If using external water, ensure the water bottle and sipper is positioned properly (Figure 1).

To open the bottle, grasp the cap and twist firmly to help separate the cap from the seal.





Ensure the feeder and water bottle hooks fully interact with the cage features.







Figure 1

Animal Care Systems | www.animalcaresystems.com | 7086 S. Revere Parkway, Centennial, CO 80112 | tel. 720.283.0177 | fax 720.283.0179

Page 1 of 1