

## Optirat GenII SOP - Rear Filter Protector for Mice



[Click for video demonstration](#)

### Introduction

The Optirat GenII filter protector guards the rear filter paper from urine when housing mice. No tools are required for installation.



**When emptying the cage, remove the filter protector assembly prior to dumping.**

### Installation

1. Remove the feeder and left rear water bottle, if necessary.
2. Insert the thin bracket onto the posts of the feeder and of the left rear water bottle; the bracket's central post must face upward (Figure Set 1).

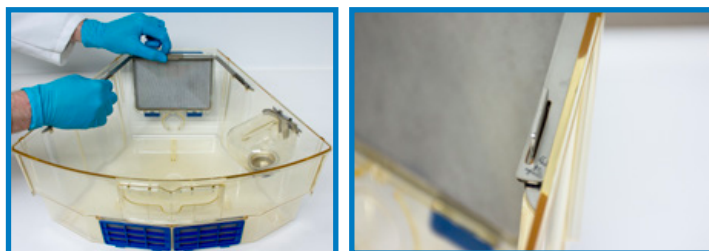


Figure Set 1

3. While stabilizing the left side of the bracket, insert the feeder onto its post (Figure Set 2).



Figure Set 2

4. Insert the protector onto its post, and ensure it's not wedged against the feeder; the guard can be shifted to the far left once installed (Figure Set 3).

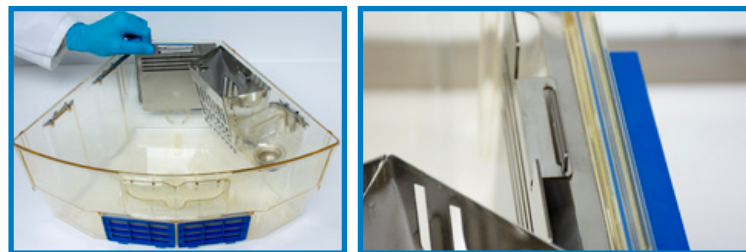


Figure Set 3



The protector is wedged against the feeder.