

Optirat GenII SOP - Upright Shelf



Video demonstration

Introductions


The Optirat GenII enables facility personnel to open a cage without bringing the cage to a biosafety cabinet.

The installation of the shelf requires a 3/16-inch Allen wrench and takes approximately five minutes.



Collar Installation

1. Remove the screws from the upright collar using a 3/16-inch Allen wrench. Insert one of the screws into the Allen wrench, and stand up the second screw for easier access for Step 3.
2. Remove the cages closest to the upright.
3. Place the two collar halves around the upright at the appropriate height, and screw them together. Stand the second screw upright so the wrench can be inserted with one hand. Ensure the screws are tight.



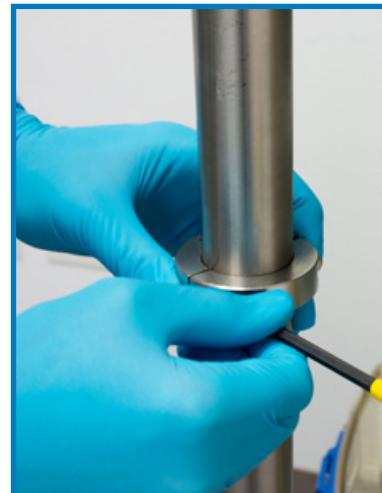


Optirat GenII SOP - Upright Shelf

Shelf Installation

1. With the shelf's hook touching or almost touching the upright, tilt the shelf to the right.
2. Push the shelf forward and pivot it so that the hook grab the shelf.
3. Lower the alignment tab into the slot into the collar.

The collar height can be adjusted with the shelf in place by loosening one or both of the screws and re-tightening.

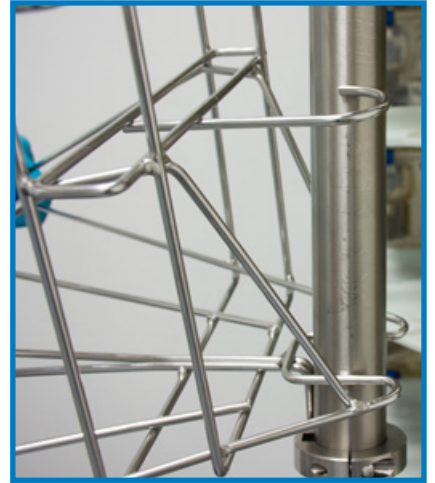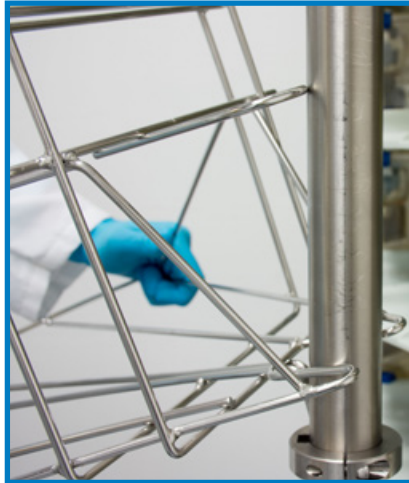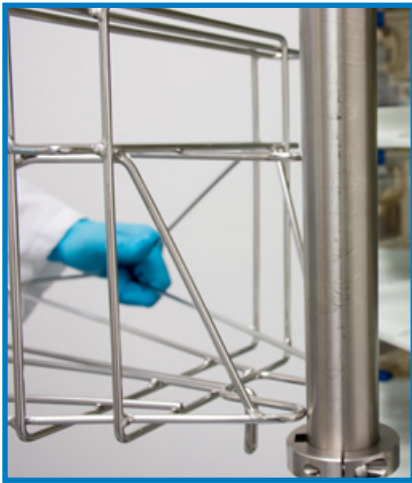

Optirat GenII SOP - Upright Shelf

Using the Shelf

1. Place an Optirat GenII cage onto the shelf.
2. Remove the cage top, and place it into the top holder narrow side first.

