


Optirat Plus SOP - Mezzanine



Video demonstration

1. Orient the mezzanine so the heads of the screws of the bottom plate face upward.

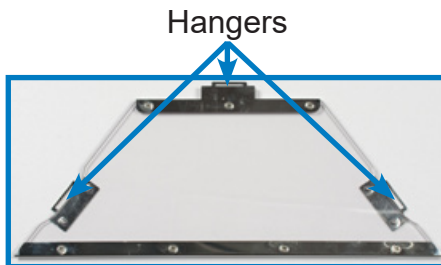

Mezzanine Sterilization Limits	
Temperature Limit	Time Limit
up to 250 F (121 C)	20 minutes



Ensure the mezzanine is installed with the proper side facing up.

2. Line up the hangers of the mezzanine with the hook features of the Optirat Plus cage base; the feeder goes beneath the mezzanine.
3. Lower the mezzanine until each hanger interacts fully with the cage's hooks (Figure 1). The hook with the feeder will accept the mezzanine hanger (Figure 2).

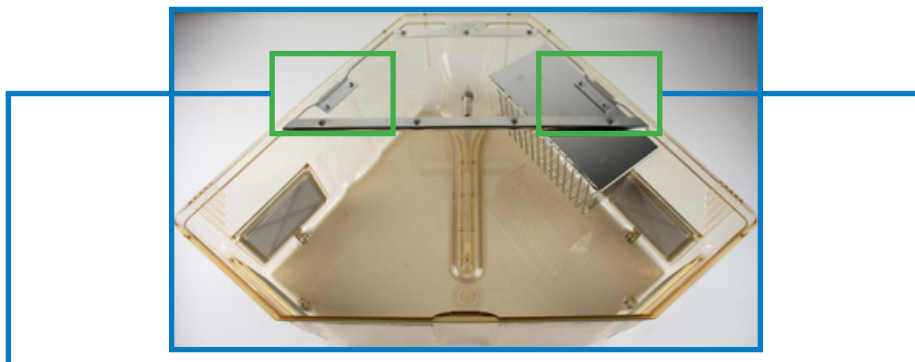


Figure 1



Figure 2